

How to use display machines

They are the computers on which we run our experiments, so we want from them the best performance. For this reason, you are not allowed to work on them, be it for writing the stimuli or analyzing data. For this you have all the other JAZZ positions. Display machines should not be online at all, if not when you need to take experiments logs. To switch from online to offline mode (which includes both JAZZ and internet), you have to click the two icons on desktop: "Enable Network", "Disable Network".

Also, the conditions under which we run our experiment must be the same, so it's crucial to check computer settings before running a session.

1. Refresh rate. Right click on desktop, -> Screen resolution -> Advanced settings -> Monitor -> Screen refresh rate
2. In the same place you can also set color settings
3. Do not accept automatic updates. Eventually report to Ayelet and she will decide.

Besides that, make sure you erase all useless data and store the good ones on DataAyelet, on JAZZ, once you're done with them.

*Lorenzo Pisoni
September 2015*